
FOOD ALLERGEN TEST STRIPS

Use these test strips for rapid detection of potential allergens in your food products. Perfect for testing solid or liquid ingredients and to prevent cross-contamination of your food items. Simple to use and results are obtain in 15 minutes. Each kit includes 25 test strips, 20 mL of dilution buffer and 80mL of extraction buffer (for gluten and soy). See accessories for a list of equipment required to prepare and manipulate your samples.

Each test is performed in 3 easy steps:

- 1- Measure 0.3mL of the liquid sample or 0.3g of the solid sample being tested (requires an optional syringe or precision balance). Solid samples need to be previously grinded to a fine powder. Add the sample to a predetermined volume of dilution buffer into a 15 mL sampling tube (see accessories). Mix the sample and the dilution buffer by shaking manually or using a vortex mixer (see accessories).
- 2- Measure 0.45mL of the extraction buffer (for gluten and soy) and add it to a small testing tube (see accessories). Add a predetermined volume of the sample mixed with the dilution buffer (0.05mL for gluten and soy or 0.5mL for casein and egg). Mix well.
- 3- Dip a test strip into the small testing tube and wait for 15 minutes. Red lines will appear to indicate either a positive or negative result.

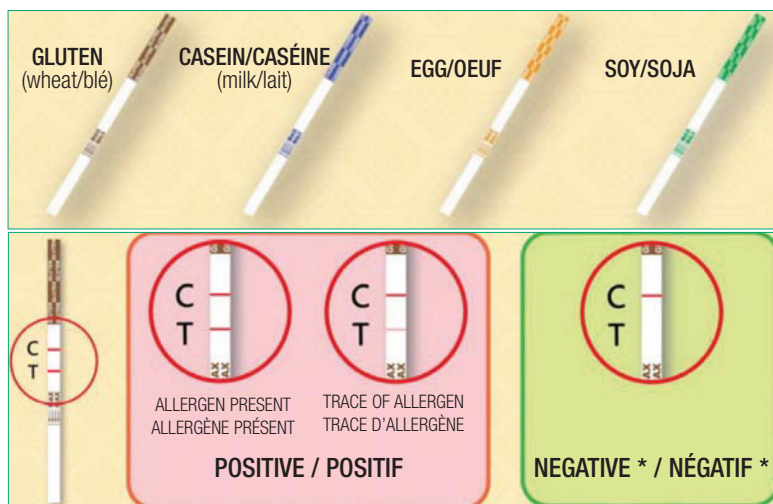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